# BILANCIA IDROSTATICA ELETTRONICA ELECTRONIC HYDROSTATIC BALANCE

GIBERTINI

## SUPERALCOMAT

laboratory  
balances

industrial  
scales

instruments  
for paint

oenological  
equipments

digital  
tensiometer

calibrated  
masses

laboratory electronic balances & oenological equipments

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**ELECTRONIC HYDROSTATIC BALANCE**  
 Fully automatic to determine Alcoholic  
 Strenght by Volume ASV, total dry  
 extract and musts evaluation.  
 Commission Regulation (EC) 2870/2000  
 and 128/2004.

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The readout of the Alcoholic Strenght by Volume is shown on a large LCD screen simultaneously with the readout of temperature, density at 20°C and other parameters.

The reading module is provided with a shower proof keyboards whose functional keys allow:

- correction of the floater volume with H<sub>2</sub>O or certified standard solutions
- zeroset and calibration of the SUPER ALCOMAT
- date, hour, minute setting and other functions
- the printing of the reading values and other parameters (printer not included)
- musts evaluation: Baumé, Babo, Brix and Oechsle degrees
- working temperature range: 10 - 30°C optimum 15 - 25°C)
- languages: Italian, German, English, French and Spanish.

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A Motorola micro-controller, specially designed for Gibertini, besides managing Super Alcomat most important functions, it is able to detect any internal problem inside the balance that could affect its accuracy, and suggest a service solution carried out by our qualified staff.

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### DENSITY

**Measurement range for the relative  
density**  
0,5 ÷ 2,25  
**Readability**  
0,00005  
**Reading repeatability**  
± 0,00005

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### Standard equipment

Power pack

2 hydrorepellent floaters, interchangeable  
in mass and volume

Container for exhaust liquids

Certified hydroalcoholic solution

### Response time

6 sec.

### Autocalibration with internal mass

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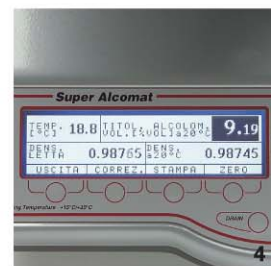
1. When the "drain" is open, flashing lamp is on 2. Automatic device to lift up and lower the floater 3. Main menu 4. Asv and density reading example 5. Device to lock the internal balance 6. Rear side: RS232 and printer outputs

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Floaters with certificate - SIT certified  
solutions 5 ÷ 60 %ASV

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### FLOATERS INTERCHANGEABLE IN MASS AND VOLUME

A particular care is devoted to the construction and calibration of the floaters.

They are all interchangeable in mass and volume to simplify their use; for example, it is possible to use two different floaters to determine the relative density of a must or a wine; in this way possible sample pollution is avoided.

In order to achieve what mentioned before, the suspension wire realised with a non-magnetic and stainless material, with a diameter less or equal to 0.20 mm. and the superficial treatment of the floater's glass were object of thorough studies to nullify the problems caused by the surface tension, by the wire's own volume, and by the electrostatic charges.

The floaters volume can, on request, be certified by competent organisations, such as the L.N.E. Paris, the I.M.G.C. Turin, the PTB Germany, etc.

### Soluzioni idroalcoliche certificate SIT dal 5% al 60% v/v - a richiesta SIT Certified hydroalcoholic solutions from 5% to 60% v/v - on request

## laboratory electronic balances & oenological equipments

customer care

info@gibertini.com

Tel. (+39) 02 3541434

ISO 9001:2000

SIT n. 94

### GIBERTINI ELETTRONICA SRL

Via Bellini, 29/31  
20026 Novate Milanese  
(Milano) - ITALY  
Tel. (+39) 02 3541434  
Fax (+39) 02 3541438  
sales@gibertini.com

www.gibertini.com